


Rees Presidio “On the Go” Web Browser

Adding New Contacts to Phone lists during alarm conditions

1. To access the Rees system, you must be on the Cleveland Clinic network (either onsite or VPN).
Type in your URL into the bar at the top of any web browser. i.e. <http://cc-clrees57> (Pharmacy).
2. Log in to the Rees system with your lab’s user name and password.

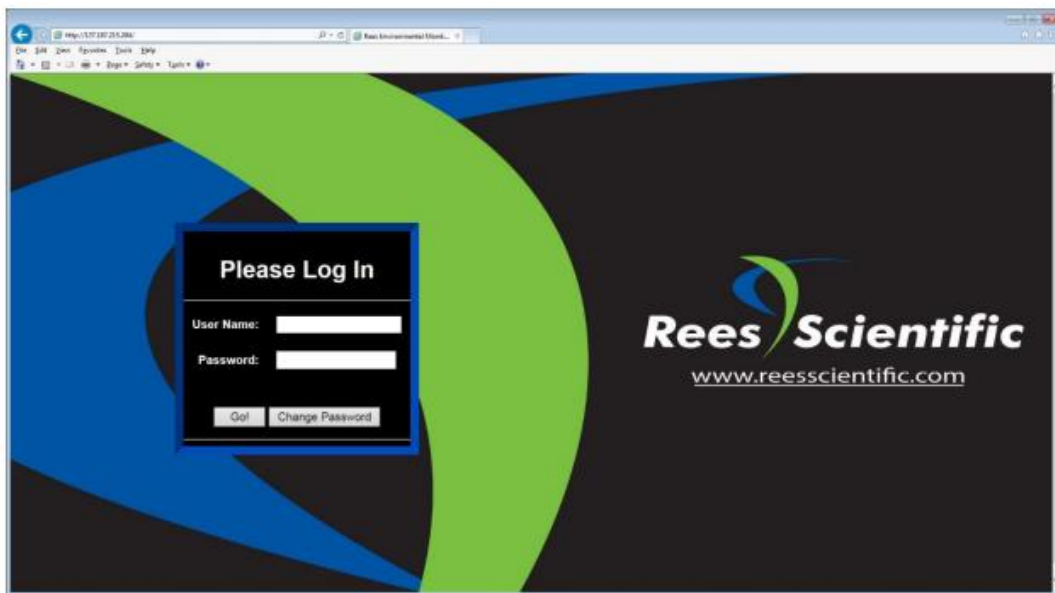

3. Select menu on the top of the status page screen




4. From the menu option select "Phone List Programming".




Adding new contacts not listed in the Phone Directory:


5. If you need to add a new emergency contact to the directory, select "Edit Directory" on the bottom of the screen.

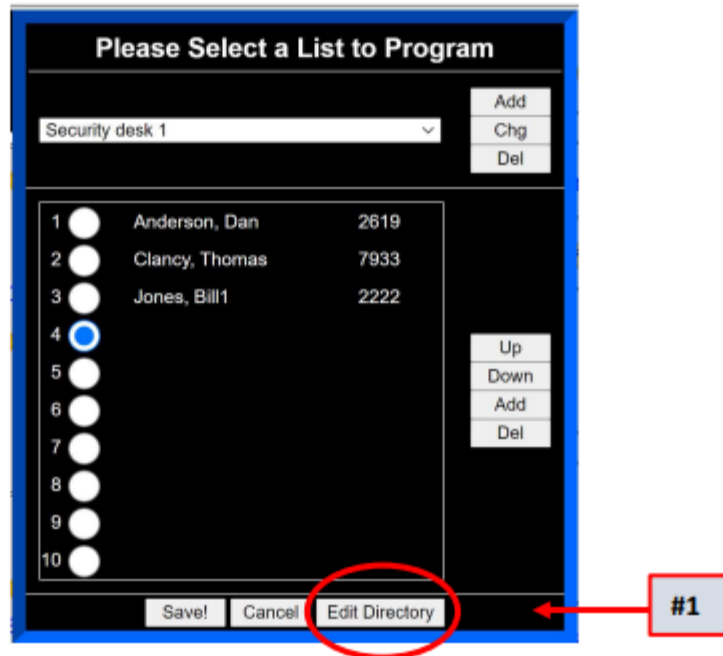

6. Select "Add" on the right-hand side of the screen.

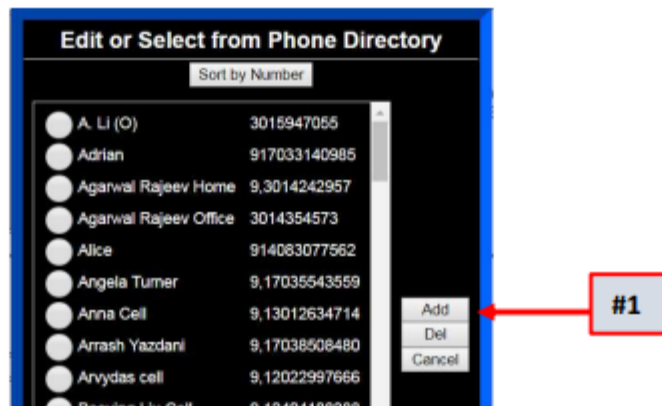

7. Enter the correct contact information in the fields and select "Save!"




Add a Phone Number

Name: Rick Saunders

Number: 609-671-3200

Save! Cancel

8. The new contact information is now in the system and can be added to any of emergency contact list from the Master directory

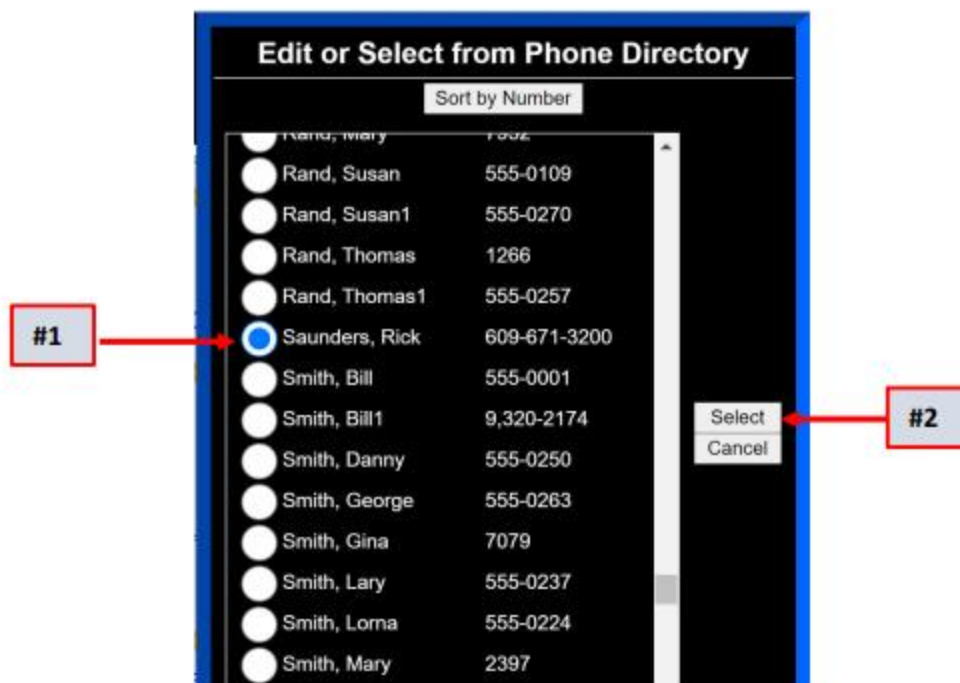

Edit or Select from Phone Directory

Sort by Number

Name	Phone
Rand, Susan	555-0109
Rand, Susan1	555-0270
Rand, Thomas	1266
Rand, Thomas1	555-0257
<input checked="" type="radio"/> Saunders, Rick	609-671-3200
Smith, Bill	555-0001
Smith, Bill1	9,320-2174
Smith, Danny	555-0250
Smith, George	555-0263
Smith, Gina	7079
Smith, Lary	555-0237
Smith, Lorna	555-0224
Smith, Mary	2397

Select
Cancel

#1

#2